



www.FreedomBreeder.com 209.664.9600 EXT. 201/202/203 Freedom Breeder · 2306 S. Daubenberger Rd. Turlock, CA 95380 USA





Product Description

This rack is made of 304 Stainless Steel, no rust ever!


RODENT 1020

Rodent 1020 racks include: 10 levels, 20 FB15 series tubs, water system, 10 attached food hoppers, base frame and casters.

These racks work perfectly for mouse holding pens, and grow out colonies; as well as African Soft Furred Rats.

Caution

Channels and backstops have sharp edges, take precautions.

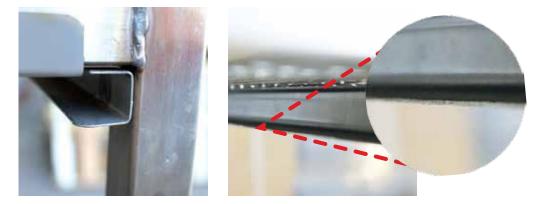

*Any comments or suggestions on this Instructions Manual please send us an email to **manuel@freedombreeder.com**



Packing List

PART #	DESCRIPTION	QTY
1	Rodent 1020 Stainless Steel Levels	10
2	FB15 Series Tubs	20
3	Water System	1
4	Attached Food Hoppers	10
5	Base Frame	1
6	4" Heavy Duty Casters (Nuts and Bolts Included)	4

*Any extra items purchased can be found on your invoice. *Packing list may vary depending on your order.

Requirements









Gloves



Two persons
*Recommended

1/2" Socket

Ratchet

Cutter





Assembly

STEP 1. Unwrapping

The Rodent 1020 is packed in a cardboard carton and plastic wrap for protection during shipment. Carefully unpack the items and remove the pallet.

Notice that the Rodent 1020 is almost assembled.

Note: If you detect any issue or damage at the moment of the delivery please contact us before you accept the shipment.



STEP 2. Parts



Continue to step 3



Assembly

STEP 3. Casters

Notice that the rack may or may not had the casters installed.


In case they have been installed, you just need to turn upside down the frame with the casters to the correct position.

In case the casters are not installed:


- 1.- Find the casters, nuts, and bolts.
- 2.- Insert the bolts into the holes in the lower frame.

3.- Using the ratchet and the 1/2" socket, tighten the caster with the bolts on each corner of the frame until the bolts are flush to the tubing of the frame.

Note: Make sure to put all the nuts and bolts included (3 per caster) and not overtighten the nuts.

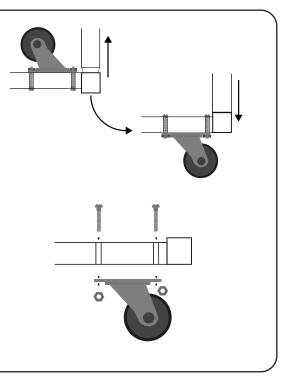

STEP 4. Stacking the levels

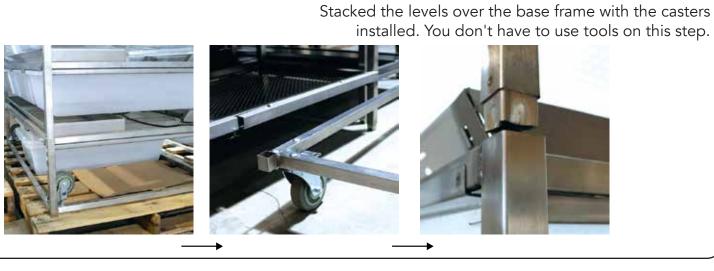





Assembly

STEP 5. Trays




STEP 6. Water system

Find the extra tube included to connect the water lines already installed.

Depending on your choice, you can use a Pressure regulator or a Gravity system.

For the **Gravity system**, put the bucket on top of your rack and let the gravity distribute the water into the water lines.

In the case of using a **Pressure regulator**, connect the water lines into the main pipe previously connected to the pressure regulator.

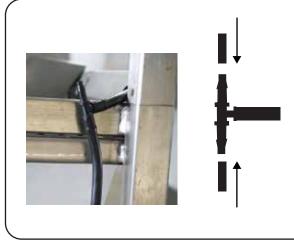


Continue to step 7







STEP 7. Water system cover (If bought)

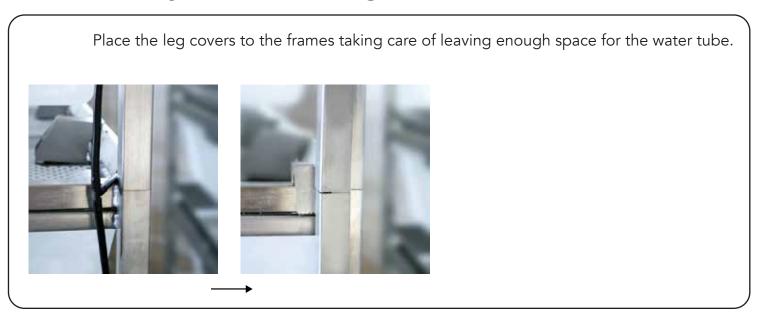


STEP 8. Water system test

- 1. Please run water through your new racking systems for a day before placing rodents or bedding in the rack.
- 2. Check for any disconnected tubing.
- 3. Adjust valves by hand to stop any leak and test again for another day. (Do not tighten with tools!)

This test is to avoid flooding your rodents.



Visit our Website, Facebook page and Instagram

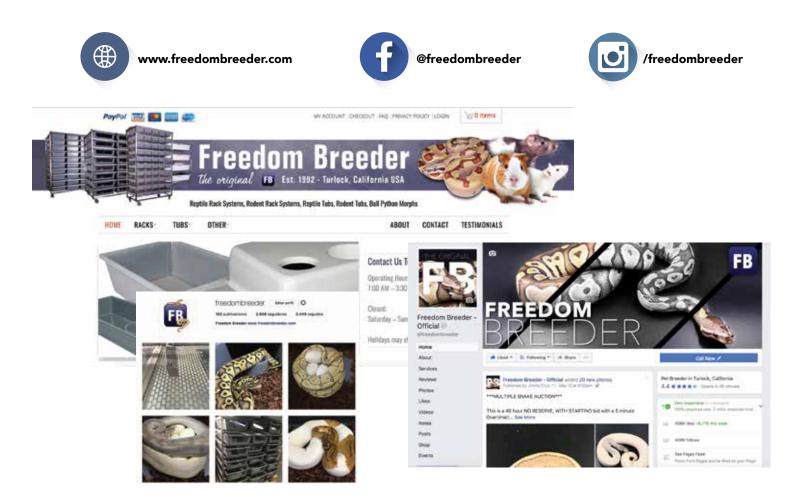

ATTENTION CUSTOMERS:

Please inform the appropriate personnel to always INSPECT YOUR SHIPMENT for damage PRIOR to signing the delivery receipt. If there is obvious damage, do not accept the shipment without NOTING IN DETAIL and taking pictures of the damage on ALL copies of the carrier's delivery receipt. Please have the freight driver SIGN ALL COPIES to acknowledge your damage notation.

-Thank You!

 *Any comments or suggestions on this Instructions Manual please send us an email to manuel@freedombreeder.com
 www.FreedomBreeder.com 209.664.9600 EXT. 201/202/203
 Freedom Breeder · 2306 S. Daubenberger Rd. Turlock, CA 95380 USA