





RODENT 99

www.FreedomBreeder.com 209.664.9600 EXT. 201/202/203 Freedom Breeder · 2306 S. Daubenberger Rd. Turlock, CA 95380 USA





Product Description


This rack is made of 304 Stainless Steel, no rust ever!

Rodent 99 racks include: 9 levels, 9 FB15 series tubs, water system, base frame and casters.

These racks work perfectly for mouse holding pens, and grow out colonies; as well as African Soft Furred Rats.

Caution

Channels and backstops have sharp edges, take precautions.

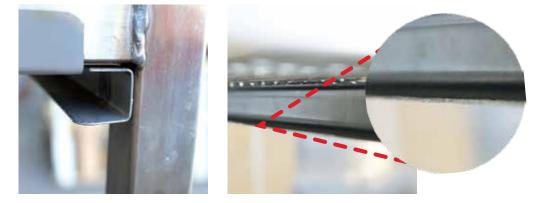

*Any comments or suggestions on this Instructions Manual please send us an email to **manuel@freedombreeder.com**

INSTRUCTION MANUAL RODENT 99 FREEDOM BREEDER



Packing List

PART #	DESCRIPTION	QTY
1	Rodent 99 Stainless Steel Levels	9
2	FB15 Semi Clear Tubs	9
3	Water System	1
4	Base Frame	1
5	4" Heavy Duty Casters (Nuts and Bolts Included)	4

*Any extra items purchased can be found on your invoice. *Packing list may vary depending on your order.

Requirements











1/2" Socket

Ratchet

Cutter

Gloves

Two persons ***Recommended**

INSTRUCTION MANUAL RODENT 99 FREEDOM BREEDER



Assembly

STEP 1. Unwrapping

The Rodent 99 is packed in a cardboard carton and plastic wrap for protection during shipment. Carefully unpack the items and remove the pallet.

Notice that the Rodent 99 is almost assembled.

Note: If you detect any issue or damage at the moment of the delivery please contact us before you accept the shipment.



STEP 2. Parts



Count all parts to make sure nothing is missing.

See Packing list for Descriptions and Quantities.

PART #	DESCRIPTION	QTY
1	Rodent 99 Stainless Steel Levels	9
2	FB15 Semi Clear Tubs	9
3	Water System	1
4	Base Frame	1
5	4" Heavy Duty Casters (Screws and Bolts Included)	4

Continue to step 3

INSTRUCTION MANUAL RODENT 99 FREEDOM BREEDER



Assembly

STEP 3. Casters

Notice that the rack may or may not had the casters installed.


In case they have been installed, you just need to turn upside down the frame with the casters to the correct position.

In case the casters are not installed:

- 1.- Find the casters, nuts, and bolts.
- 2.- Insert the bolts into the holes in the lower frame.


3.- Using the ratchet and the 1/2" socket, tighten the caster with the bolts on each corner of the frame until the bolts are flush to the tubing of the frame.

Note: Make sure to put all the nuts and bolts included (3 per caster) and not overtighten the nuts.

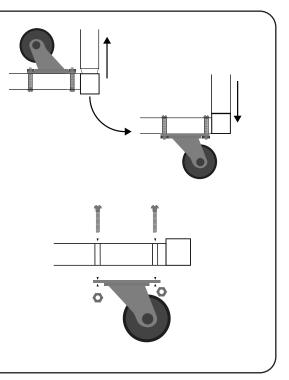

STEP 4. Stacking the levels

Stacked the levels over the base frame with the casters installed. You don't have to use tools on this step.

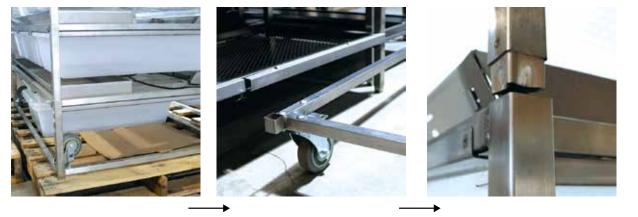






Assembly

STEP 5. Trays




STEP 6. Water system

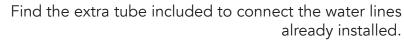

Depending on your choice, you can use a Pressure regulator or a Gravity system.

For the **Gravity system**, put the bucket on top of your rack and let the gravity distribute the water into the water lines.

In the case of using a **Pressure regulator**, connect the water lines into the main pipe previously connected to the pressure regulator.

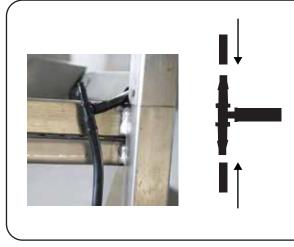


Continue to step 7







STEP 7. Water system cover (If bought)

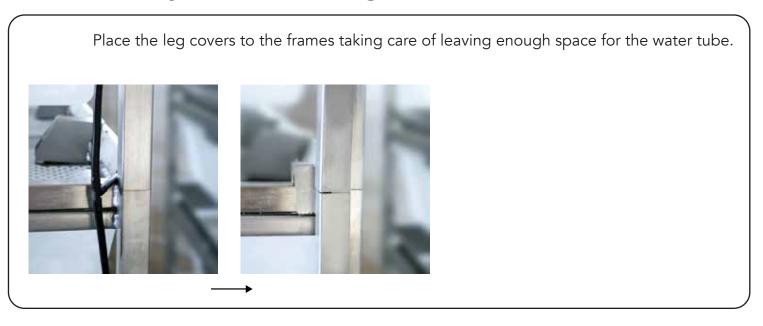


STEP 8. Water system test

- 1. Please run water through your new racking systems for a day before placing rodents or bedding in the rack.
- 2. Check for any disconnected tubing.
- 3. Adjust valves by hand to stop any leak and test again for another day. (Do not tighten with tools!)

This test is to avoid flooding your rodents.



Visit our Website, Facebook page and Instagram

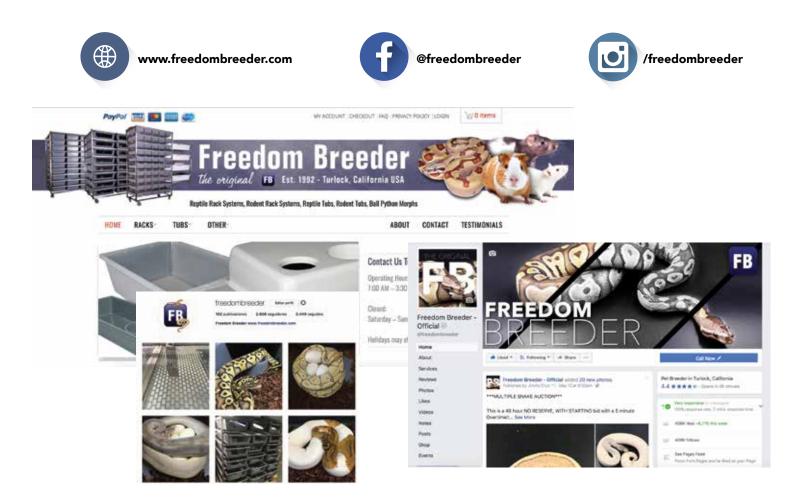

ATTENTION CUSTOMERS:

Please inform the appropriate personnel to always INSPECT YOUR SHIPMENT for damage PRIOR to signing the delivery receipt. If there is obvious damage, do not accept the shipment without NOTING IN DETAIL and taking pictures of the damage on ALL copies of the carrier's delivery receipt. Please have the freight driver SIGN ALL COPIES to acknowledge your damage notation.

-Thank You!

 *Any comments or suggestions on this Instructions Manual please send us an email to manuel@freedombreeder.com
www.FreedomBreeder.com 209.664.9600 EXT. 201/202/203
Freedom Breeder · 2306 S. Daubenberger Rd. Turlock, CA 95380 USA